


MAY 2020

## *High Pressure Reactor Improves Development of Novel Catalysts*

### **The Chianese Organometallic Chemistry group**

at Colgate University (Hamilton, NY, USA) are using a custom designed Multicell Parallel High-Pressure Reactor from Asynt to conduct reproducible hydrogenation experiments in parallel, to rapidly determine the kinetic properties of their novel catalysts.

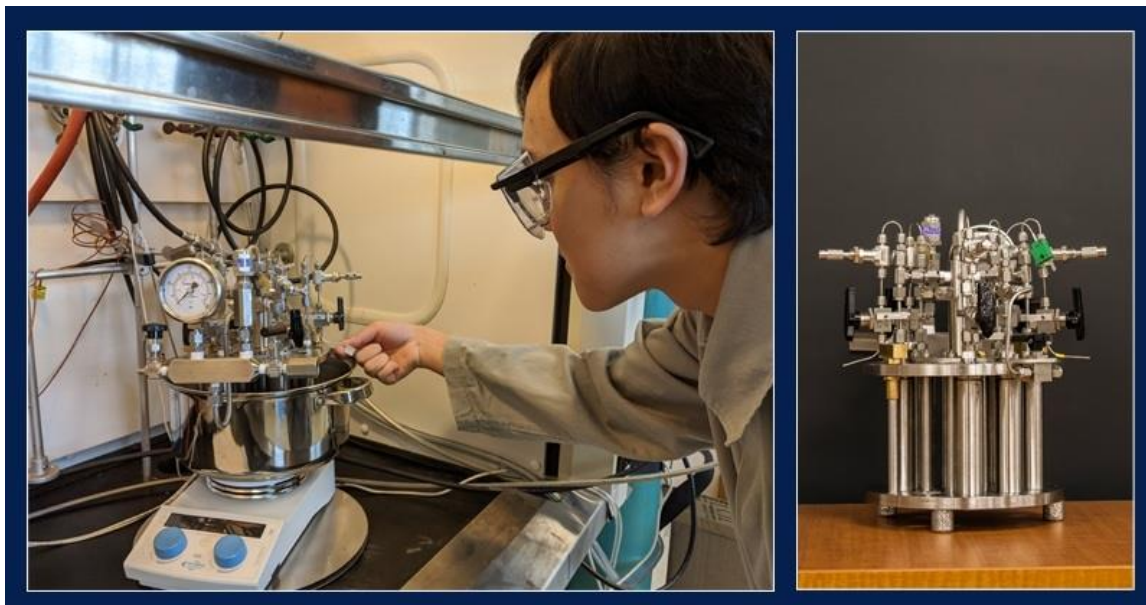

### **The research activities**

of the Chianese group focuses on the development of new transition metal complexes for catalysing the hydrogenation of polar compounds. To accelerate this process the group wanted a high-pressure reactor capable of conducting many hydrogenation reactions at once, with rigorous control of the initial concentrations and reaction temperature. They also needed the ability to take and analyse reaction aliquots over the time course of the reaction, which could then be used to for kinetic profiling.

### **Associate Professor Anthony Chianese,**

head of the group, commented “By working with the technical team at Asynt we were able to specify a custom version of their Multicell parallel high-pressure reactor that exactly fitted our needs. Before we had the Multicell we were running several test tube reactions within a single batch reactor, meaning that we could only conduct analysis at the end and would often experience contamination between samples. The Multicell now allows us to run many reactions in parallel

#### **Asynt Ltd**

Unit 29 Hall Barn Road Industrial Estate Isleham Cambridgeshire United Kingdom CB7 5RJ  
T: +44 (0)1638 781709 F:+44(0)1638 781706 sales@asynt.com www.asynt.com

Registered office: Eldo House, Kempson Way, Bury St Edmunds, Suffolk, IP32 7AR Registration No: 5160407  
VAT No: GB 838 5592 82



with no detectable crossover between the headspace of different cells. Importantly, because we can now analyse aliquots while the reactions are running, it is possible to rapidly determine reaction rates for any given catalytic hydrogenation. Often our organometallic catalysts are highly air sensitive, so it is great to be able to assemble our Multicell reactor inside a glovebox then bring it out to run the reactions. We also really like the fact that the whole unit is user-serviceable with all of the parts, aside from the cells themselves, easily replaceable if there is an issue.”

#### **When asked if he would recommend the system**

to other researchers, Professor Chianese responded “Yes. I think most researchers only analyze hydrogenation reactions at the endpoint because they don’t have a way to take aliquots. With the Asynt Multicell you can study hydrogenation reactions much more quickly, efficiently and in more detail than was previously possible.”

#### **The Multicell high pressure reactor**

is designed for undertaking experiments including carbonylation, hydrogenation, catalysis and supercritical fluid reactions, as well as air sensitive material studies. Used in conjunction with a hotplate stirrer the standard 316 stainless steel Multicell reactor offers users the ability to undertake 10 x 30 mL reactions in parallel at pressures of up to 50 bar and temperatures up to 200 °C. To increase the versatility of this high-performance system, Asynt can work with the end user on specific customisation to enhance the capabilities of the Multicell towards certain applications. Options include the ability to make additions/take samples under pressure, increased operating conditions up to 350 °C and 200 bar, a 4-position system offering reaction volumes of up to 150 mL, sub-ambient reactor cooling, and the ability to have different temperatures and pressures in each reactor cell.

#### **For further information on**

the Multicell Parallel High Pressure Reactor please visit <https://www.asynt.com/product/multicell-parallel-high-pressure/> or contact Asynt on +44-1638-781709 / [enquiries@asynt.com](mailto:enquiries@asynt.com). For further information on the work of the Chianese Group at Colgate University please visit <https://www.colgate.edu/about/directory/achianese>

#### **Asynt Ltd**

Unit 29 Hall Barn Road Industrial Estate Isleham Cambridgeshire United Kingdom CB7 5RJ  
T: +44 (0)1638 781709 F:+44(0)1638 781706 sales@asynt.com www.asynt.com

Registered office: Eldo House, Kempson Way, Bury St Edmunds, Suffolk, IP32 7AR Registration No: 5160407  
VAT No: GB 838 5592 82



## **Asynt**

is a leading supplier of affordable products, consumables and services for chemists in industry and academia. With staff of trained chemists - Asynt can draw upon this in-depth applications knowledge to provide a high level of customer support for its DrySyn Heating Blocks, Controlled Lab Reactors, Synthesis Tools, Evaporators, Circulators, Temperature Control Systems, Vacuum Pumps and Lab Safety Equipment.

## **Worldwide HQ**

### **Asynt Ltd**

Unit 29, Hall Barn Industrial Estate  
Isleham  
Cambridgeshire CB7 5RJ  
UK

tel: +44-1638-781709

email: [enquiries@asynt.com](mailto:enquiries@asynt.com)

web: [www.asynt.com](http://www.asynt.com)

### **Asynt Ltd**

Unit 29 Hall Barn Road Industrial Estate Isleham Cambridgeshire United Kingdom CB7 5RJ  
T: +44 (0)1638 781709 F:+44(0)1638 781706 sales@asynt.com www.asynt.com

Registered office: Eldo House, Kempson Way, Bury St Edmunds, Suffolk, IP32 7AR Registration No: 5160407  
VAT No: GB 838 5592 82